


Kirchhoff's laws

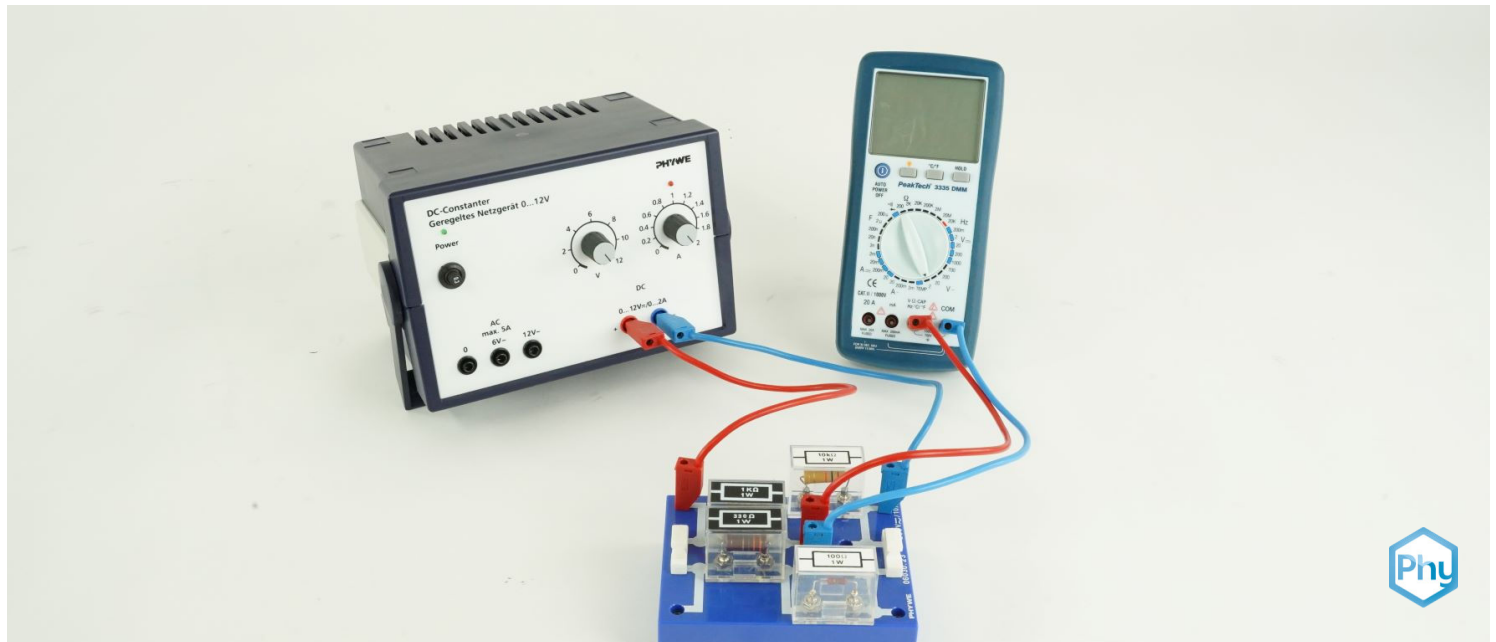

Physics

Electricity & Magnetism

Electric current & its effects

Physics

Electricity & Magnetism

Simple circuits, resistors & capacitors

Applied Science

Engineering

Electrical Engineering

Properties of Electrical Circuits



Difficulty level

hard



Group size

2



Preparation time

45+ minutes



Execution time


45+ minutes

This content can also be found online at:



<http://localhost:1337/c/621607dc83e26600034d676d>

PHYWE

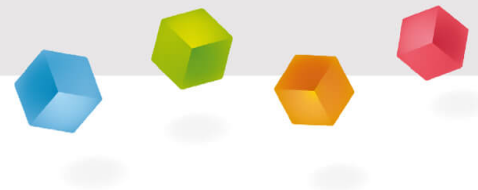

General information

Application

PHYWE



Fig. 1: Experimental set-up

The Kirchhoff's laws are to equalities which deal with the current in a node and voltage around a loop. These two laws and the equations for individual components are the basis for analyzing circuits.

Other information (1/2)

PHYWE

**Prior****knowledge****Main****principle**

The prior knowledge for this experiment is found in the Theory section.

First Kirchhoff's laws are verified by measuring current, voltage and resistance in series and parallel circuits. The Wheatstone bridge circuit is used to determine unknown resistances more precisely.

Other information (2/2)

PHYWE

**Learning****objective****Tasks**

The goal of this experiment is to analyze circuits using the two Kirchhoff's laws.

1. Task: Verify Kirchhoff's laws by measuring current and voltage for series and parallel connected resistors for each resistor as well as the total values. From these measurements calculate the partial and total resistances.

2. Task: Determine unknown resistances by the use of the Wheatstone bridge circuit.

Theory (1/3)

PHYWE

Task 1: With branched circuits, in the steady-state condition, Kirchhoff's 1st law applies at every junction point:

$$\sum_k I_k = 0 \quad (1)$$

where I_k are the currents leading to or from the junction point. This means, that in every junction point, the charge is conserved. It is customary to take I_k as negative if the corresponding current in the k -th conductor is flowing away from the junction point. For every closed loop C in a network of linear conductors, in the steady-state condition, Kirchhoff's 2nd law applies:

$$\sum_k U_k = 0 \quad (2)$$

where U_k the voltage in the k -th conductor. This is a special case of the induction law as it applies only for constant magnetic flows. More precisely, it is a conclusion of the 1st and 3rd of Maxwell's equations. It means that, in a closed loop, the electrical energy is conserved.

Theory (2/3)

PHYWE

From these laws follow some conclusions how current, resistance and voltage behave in parallel and series circuits:

1. series connection of several conductors

$I_n = I_m$: Through every conductor flows the same current.

$U_{tot} = \sum_n U_n$ The voltages across the individual conductors sum up to the total voltage across the circuit.

$R_{tot} = \sum_n R_n$ The resistances across the individual conductors sum up to the total resistance across the circuit.

From this follows: $\frac{U_{tot}}{R_{tot}} = \frac{U_n}{R_n}$ as well as $\frac{R_n}{R_{tot}} = \frac{U_n}{U_{tot}}$ (3)

where $I = \frac{U}{R}$ is Ohm's law.

Theory (3/3)

PHYWE

Task 2: In principal an unknown resistance can be determined by measuring current and voltage across the resistance. The finite intrinsic resistances of the instruments would introduce significant errors. To avoid such errors the measurement has to be done current-free. In a Wheatstone bridge circuit the unknown resistance R_x is connected to three resistances (see Fig. 4) of which at least one is variable.

In this experiment R_3 is adjusted in such a manner that no current is flowing through the instrument G (alignment of the bridge) which means the voltage across G vanishes as well. In this case the voltages across R_x and R_1 are equal as well as across R_3 and R_2 . Also, as no current is flowing through G , the same current is flowing through R_3 and R_x on the one hand (denoted as I_1) and through R_1 and R_2 on the other hand (I_2). This results into the following equations:

$$I_1 R_x = I_2 R_1 \text{ and } I_1 R_3 = I_2 R_2$$

Theory (4/4)

PHYWE

Division of these relations yields eq. (4) which computes the unknown resistance R_x .

$$R_x = R_3 \cdot \frac{R_1}{R_2} \quad (4)$$

The voltage at the source is unimportant and may even be time-dependent.

Note: In a customary Wheatstone bridge, R_1 and R_2 are two parts of a wire separated by a sliding contact (similar to a potentiometer). So R_1 and R_2 are both variable where the sum is the total resistance of the wire. However, in this experiment R_1 and R_2 are constant and R_3 is changed.

Equipment

Position	Material	Item No.	Quantity
1	PHYWE Power supply, 230 V, DC: 0...12 V, 2 A / AC: 6 V, 12 V, 5 A	13506-93	1
2	Digital multimeter, 600V AC/DC, 10A AC/DC, 20 MΩ, 200 μF, 20 kHz, -20°C... 760°C	07122-00	1
3	Resistor 100 Ohm, 1W, G1	39104-63	1
4	Resistor 220 Ohm, 1W, G1	39104-64	1
5	Resistor 330 Ohm, 1W, G1	39104-13	1
6	Resistor 470 Ohm, 1W, G1	39104-15	1
7	Resistor 1 kOhm, 1W, G1	39104-19	2
8	Resistor 2.2 kOhm, 1W, G1	39104-23	1
9	Resistor 3.3 kOhm, 1W, G1	39104-25	1
10	Resistor 4.7 kOhm, 1W, G1	39104-27	1
11	Resistor 10 kOhm, 1W, G1	39104-30	1
12	Connection box	06000-00	1
13	Short-circuit plug, white	06027-06	3
14	Connecting cord, 32 A, 250 mm, red	07360-01	2
15	Connecting cord, 32 A, 250 mm, blue	07360-04	2

PHYWE



Setup and Procedure

Setup (1/2)

PHYWE

Task 1: The circuit is set up as shown in Fig. 2. The digital multimeter has to be used as Voltmeter and Ampèremeter alternatingly. Try different resistors in order to verify Kirchhoff's laws.



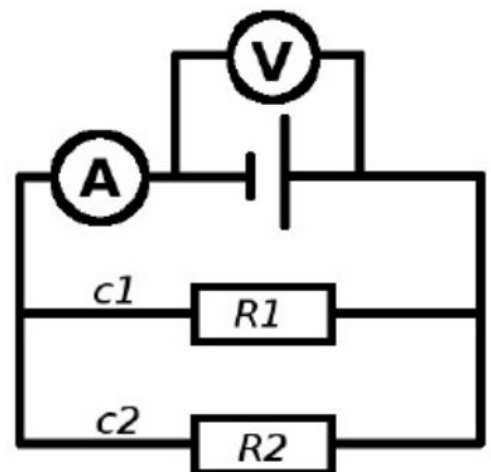


Fig. 2: Schematic circuit for Task 1.

Setup (2/2)

PHYWE

Task 2: The resistors that have to be determined need to be disguised. You can use non-transparent tape to cover the resistors' cases. Set up the experiment according to the figures on the right. The digital multimeter is connected in series with and in order to measure the current. Choose the lowest possible measuring range and set the measuring mode to direct current (DC). Be careful to plug the connection cords into the correct sockets of the multimeter. If actually using a Wheatstone bridge, R_1 and R_2 are variable with $R_1 + R_2 = \text{const.}$

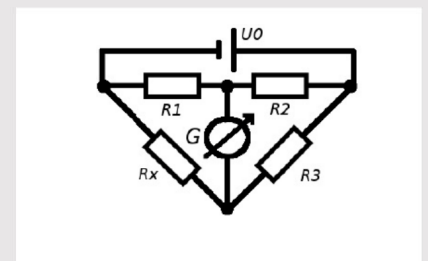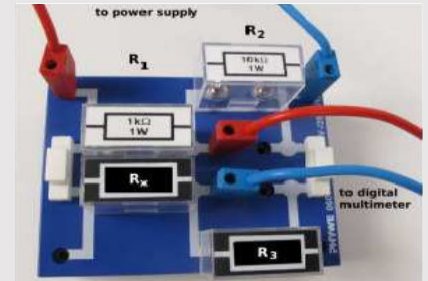


Procedure

PHYWE

Task 1: Before switching on the power supply, make sure that both adjustments of current and voltage are tuned down to zero. After switching on the power supply first tune the current until the green LED goes out. Then carefully tune up the voltage to a maximum of 3V. Measure the current in the unbranched part of the circuit. Replace $c1$ with the multimeter and measure the partial current I_1 . Continue with the measurement of I_2 at $c2$. Measure the applied voltage directly at the source. Afterwards, insert a third resistor R_3 in series with R_1 . Measure the partial voltages U_1 and U_3 across R_1 and R_3 respectively. Note: To perform the measurements you will have to build the circuit for each measurement individually.

Task 2: Before switching on the power supply, make sure that both adjustments of current and voltage are tuned down to zero. After switching on the power supply first tune the current until the green LED goes out. Then carefully tune up the voltage to a maximum of 1V. If the green LED lights up again you have to adjust the current. Keep an eye on the digital multimeter measuring the current in the circuit and keep the current well below 1A. In order to determine the unknown resistance change R_3 until the current through G vanishes. You may try single resistors as well several resistors connected in series to vary R_3 . Note down the combinations at which the current becomes zero or at least reaches a minimum.

PHYWE

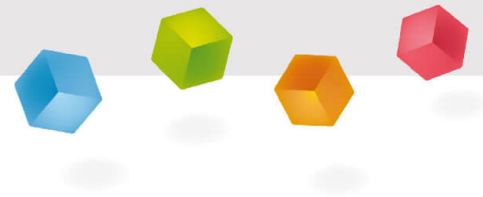

Evaluation

Task 1

PHYWE

In the following the evaluation of the obtained values is described with the help of example values. Your results may vary from those presented here.

In order to verify Kirchhoff's laws compare the measured values with the theoretical values obtained from eqs. (1) to (3). To some degree, deviations from the computed results have to be expected, e.g. the connection cords have a non-zero resistance and therefore contribute to the total resistance of the circuit.

Task 2

PHYWE

Having found the appropriate resistances for R_3 that will minimize the current through G use eq. (4) to calculate the unknown resistances. Tab 1 shows an example of combinations of R_3 and R_x for the setup of this experiment.

	R_3 [$k\Omega$]	R_x [$k\Omega$]
X1	1.1	0.1
X2	2.2	0.22
X3	3.3	0.33
X4	4.7	0.47

Tab. 2: Resistances determined via the Wheatstone bridge circuit with $R_1 = 1k\Omega$ und $R_2 = 10k\Omega$

Instead of single resistors you can use several potentiometer with different ranges connected in series or a decade resistance box (06194-10) as the third resistance R_3 .